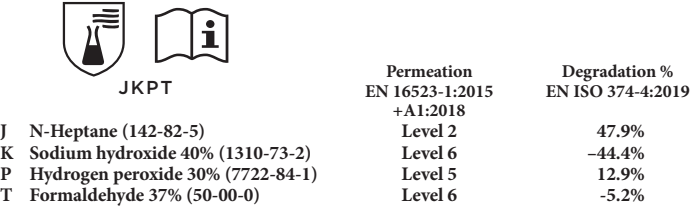
## EN ISO 374-1:2016+A1:2018/Type B

|                  |                   |                   |
|------------------|-------------------|-------------------|
| Level 1 > 10 min | Level 3 > 60 min  | Level 5 > 240 min |
| Level 2 > 30 min | Level 4 > 120 min | Level 6 > 480 min |

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[EN]: Donning: Thoroughly wash hands. Select the appropriately sized gloves. Hold with one hand and insert the other. Pull glove cuff towards wrist to cover as much skin as possible and secure glove. Check to make sure there are no holes or tears. Doffing: Grasp the outside edge of the glove near the wrist. Peel the remaining glove off from the inside, creating a "bag" containing both gloves. Peel the glove away from the hand, turning it inside out.

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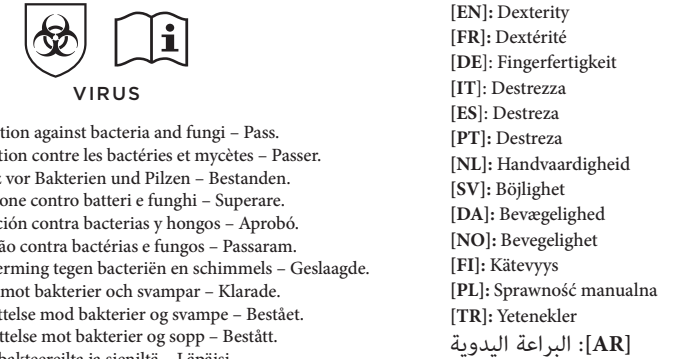
EN ISO 21420:2020

|         |       |
|---------|-------|
| Level 1 | 11mm  |
| Level 2 | 9,5mm |
| Level 3 | 8mm   |
| Level 4 | 6,5mm |
| Level 5 | 5mm   |

| Performance level | AQL   | Inspection levels |
|-------------------|-------|-------------------|
| Level 3           | <0,65 | G1                |
| Level 2           | <1,5  | G1                |
| Level 1           | <4,0  | S4                |

ChemRest Series EBT™ by Showa: The enhanced EBT Technology (EBT) are engineered for applications in laboratories performing long-term landfill leachate testing. Showa's gloves with EBT achieved 82.0% reduction in 1.9% biodegradation over the same period.

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Gloves have a minimum shelf life of 5 years.  
Gloves provide protection from chemical hazards shown.  
Potential allergens: carbamate

The penetration resistance has been assessed under laboratory conditions to the tested specimen.

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It is recommended to check that the gloves are suitable for the intended conditions at the workplace may differ from the type test depending on abrasion and degradation. When used, protective gloves may provide protection against dangerous chemical due to changes in physical properties. Movement of the gloves may cause degradation caused by the chemical contact etc. may reduce the actual protection.

Before usage, inspect the gloves for any defects or imperfections. All gloves should be stored unopened at room temperature and protected from heat, humidity, sunlight, ozone, pests and sharp objects. It is recommended that all Showa products be used in accordance with the instructions on the label.

venleri seçin. Bir elinizle tutun ve First in, First Out stock rotation. Products with or without an expiration date. Do not use expired products. Discard the customer from ensuring product viability upon receipt and do not use a product for any particular use.

veya yırtık olmadığından emin olun. Discard used gloves in compliance with local regulations. Do not wear gloves that are damaged or worn. Discard gloves that are damaged or worn. Discard gloves that are damaged or worn.

divenin geri kalanını soyarak a risk of entanglement by moving parts of a machine.

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